I 12 INCONVENIENTI E RIMEDI

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GB 12 TROUBLESHOOTING

The maintenance operations that ANGELANTONI INDUSTRIE authorizes are those indicated in the chapter entitled "Regular maintenance". The information given in the column "SOLUTION" in the tables below does not authorize any operations if these compromise safety; the information is given to help any specialized technicians find the fault.

PROBLEM	PROBABLE REASON	SOLUTION
The machine does not start.	<ul style="list-style-type: none"> There is no power supply from the main. The main switch is pointing to "OFF". The ON/OFF switch is pointing to "OFF". A mistake in the "RUN" programming. Alarm signal on the display. Some other reason not mentioned above. 	<ul style="list-style-type: none"> Check and proceed accordingly. Turn on the main switch. Turn on the ON/OFF switch. Consult KEYKRATOS™ PLUS sw and/or the WINKRATOS™ sw handbooks and proceed accordingly. Check the description of the type of alarm on the LCD display and eliminate the cause. Call the technical assistance service.
The equipment does not reach the set temperature.	<ul style="list-style-type: none"> Error in setting of setpoint. There is too much material in the test chamber. The test material prevents correct air circulation. The test material dissipates too much heat. The door is not shut correctly. 	<ul style="list-style-type: none"> Check and modify. Reduce the quantity. Place the material so that air can circulate through the intake and outlet openings. Check and reduce the load that dissipates the heat according to the technical specifications of the chamber. Check that nothing is preventing it from closing.

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